3326KR  
Lot 002~

Printed February 04, 2020  
Version 1.1

For Research Use Only. Not for use in diagnostic procedures.

## **DDDDK-tagged Protein PURIFICATION CARTRIDGE**

(Clone FLA-1GS)

CODE No. 3326KR

*PURIFICATION to maintain protein activity  
from eukaryote cell lysate and culture supernatant*

## **Product Description**

Recombinant tagged proteins are widely used for protein purification and protein characterization. There are many purification methods and kits available. DDDDK epitope tag (DYKDDDDK) has been widely used as a multi-purpose tag for protein purification.

MBL's DDDDK-tagged Protein PURIFICATION CARTRIDGE is ready-to-use-cartridge for the isolation of DDDDK-tagged protein from cell culture supernatants and cell lysate under neutral pH condition. The cartridge can be used directly on automated ÄKTA FPLC Systems.

The elution of DDDDK-tagged proteins from the cartridge is achieved by the addition of the DDDDK-tag peptide (DYKDDDDK). As the DDDDK-tag peptide competes with DDDDK-tagged proteins on the cartridge, the purified DDDDK-tagged proteins maintain function and retain protein activity.

## **Components**

	Quantity
	CODE No. 3326KR
<b><u>DDDDK-tagged Protein PURIFICATION CARTRIDGE</u></b>	<b><u>1 mL × 1</u></b>
• DDDDK-tag cartridge (1/16 inch female end) × 1, Stop plugs (1/16 inch male) × 2	

Elution Peptide (DYKDDDDK) is NOT included.

It is available as CODE No. 3325-205 (1 mg × 5 vials).

## **Storage**

Store for up to 1 year from date of receipt at 2-8°C. Do not freeze.

## Cartridge Specifications

<b>Specifications</b>	
Column volume	1 mL
Cartridge body material	Polypropylene
Connector size	1/16 inch
Matrix	Cross-linked agarose
Ligand	Anti-DDDDK-tag mAb (clone: FLA-1GS)
Binding capacity	Approx. 2 mg recombinant DDDDK-tagged Protein/mL gel
Maximum pressure	0.3 MPa
Recommended flow rate	0.5-1 mL/min
Storage buffer	0.1% ProClin 150/PBS

## Material Preparation

Prepare the following reagents before affinity purification.

1. Lysis buffer : Suitable Lysis buffer varies with a kind of the DDDDK-tagged protein  
(See **Additional Information**).  
[Example of buffer constitution]  
10-50 mM Tris-HCl (pH 7.5)  
100-300 mM NaCl  
1% NP-40 or Triton X-100  
If necessary add Protease Inhibitor Cocktail  
(e.g. SIGMA: code P8340, PIERCE: code 78415)
2. Washing buffer : Suitable Washing buffer varies with a kind of the DDDDK-tagged protein  
(See **Additional Information**).  
[Example of buffer constitution]  
10-50 mM Tris-HCl (pH 7.5) or HEPES-KOH (pH 7.5)  
300-500 mM NaCl  
To prevent non-specific protein binding to the gel,  
salt concentration should be 300-500 mM NaCl.
3. Elution buffer : 0.1 mg/mL DDDDK-tag peptide in PBS (or Washing buffer).
4. Regeneration buffer : 0.17 M Glycine-HCl (pH 2.3)
5. Column storage buffer : 0.1% ProClin 150/PBS
6. A suitable liquid chromatography system with 1/16 inch tubing (e.g. GE Healthcare: ÄKTA system)

Note: For best results, filter all the buffers through 0.45 µm filter and degas before use.

## **Protocols**

### A. Sample preparation

The cartridge is optimized under only native conditions. It is not recommended for use under denaturing conditions or for purification of aggregated, unstable, and insoluble protein (e.g. inclusion bodies). Proteins solubilized with such as 6 M Guanidine-HCl or 8 M Urea cannot be purified using this cartridge (see **Additional Information**).

Cellular debris and particulate matter must be removed prior to purification. The protein extract should be centrifuged (10,000-20,000 × g for 15 min) and filtered with a 0.45 µm filter to remove any remaining cells and particulates.

Highly viscous samples containing chromosomal DNA or RNA should be sonicated or treated with nuclease to reduce viscosity.

### B. Purification

1. Fill the pump A and B in liquid chromatography system with Washing buffer and Elution buffer.
2. Remove one end of the stop plug from the cartridge, and start pumping Washing buffer at a flow rate of 0.5 mL/min until a few drops fill in the top of the cartridge. Connect the cartridge to the pump side tubing.
3. Immediately remove the other end of the stop plug from the cartridge, connect the cartridge to the UV monitor side tubing.
4. Equilibrate the cartridge with 5-10 bed volumes of the Washing buffer at a flow rate of 1 mL/min.
5. Apply the sample to the cartridge at a flow rate of 0.5-1 mL/min. Collect fractions.
6. Wash the cartridge with 20 bed volumes of the Washing buffer until the absorbance approaches baseline at a flow rate of 1 mL/min.
7. Elute with 8 bed volumes of Elution buffer and collect 0.5 mL fractions.

### C. Regeneration and storage

1. Wash the cartridge with 10 bed volumes of Regeneration buffer.
2. Immediately wash the column with 10 bed volumes of Column storage buffer.
3. Screw the stop plug tightly, store the cartridge at 2-8°C.



Note: This cartridge may be used multiple times, depending on the usage conditions.

## **Related Products:**

Please visit our website at <https://ruo.mbl.co.jp/>.

## Example of Purification Result

### Purification of N-terminus DDDDK-tagged $\beta$ -Galactosidase

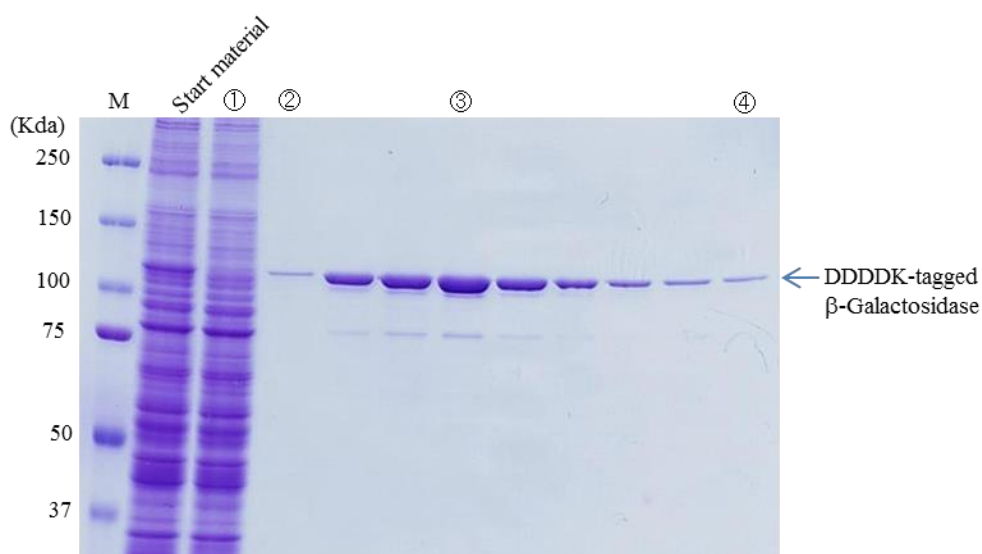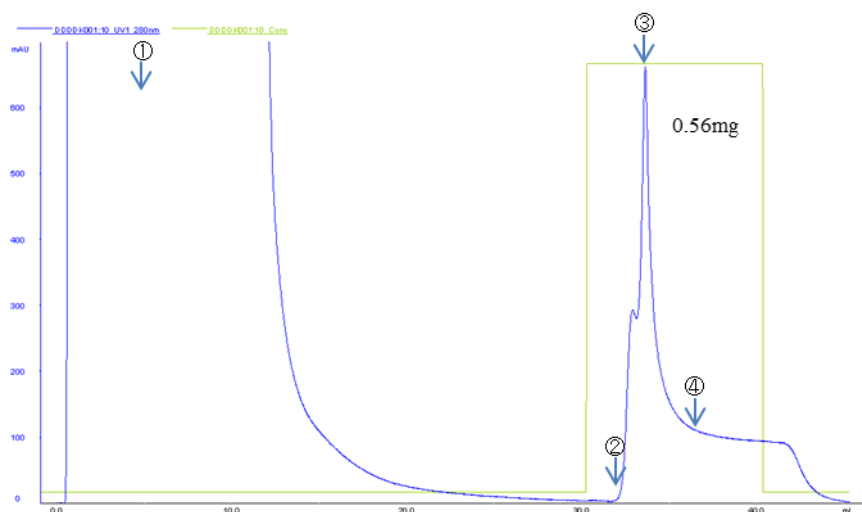Cartridge	: DDDDK-tagged Protein PURIFICATION CARTRIDGE
Sample	: DDDDK-tagged $\beta$ -Galactosidase/293T transient transfectant ( $1 \times 10^8$ cells)
Cell lysis buffer	: 10 mM Tris-HCl, 150 mM NaCl, 1% NP-40 (pH 7.5)
Wash buffer	: 10 mM Tris-HCl, 300 mM NaCl (pH 7.5)
Elution buffer	: 0.1 mg/ml DYKDDDDK peptide in PBS
Flow rate	: 1.0 mL/min (0.5 mL/min for elution)
Chromatography system	: ÄKTAexplorer 10S (GE Healthcare)

## **Additional Information**

Several reagents were examined whether or not they were suitable for use with the DDDDK-tagged Protein PURIFICATION CARTRIDGE. For example, RIPA buffer could be used for preparation of cell lysate. The results are listed below.

### **Chaotropic agents**

<b>Urea</b>	<b>1 M</b>	<b>Yes</b>
<b>Guanidine-HCl</b>	<b>1 M</b>	<b>No</b>

### **Reducing agents**

<b>DTT</b>	<b>10 mM</b>	<b>Yes</b>
<b>2-Mercaptoethanol</b>	<b>10 mM</b>	<b>Yes</b>

### **Surfactants**

<b>Nonionic</b>	<b>Tween-20</b>	<b>5%</b>	<b>Yes</b>
	<b>TritonX-100</b>	<b>5%</b>	<b>Yes</b>
	<b>NP40</b>	<b>1%</b>	<b>Yes</b>
	<b>Digitonin</b>	<b>1%</b>	<b>Yes</b>
	<b>n-Octyl-<math>\beta</math>-D-gulcoside</b>	<b>1%</b>	<b>Yes</b>
<b>Zwitterionic</b>	<b>CHAPS</b>	<b>1%</b>	<b>Yes</b>
	<b>CHAPSO</b>	<b>1%</b>	<b>Yes</b>
<b>Anionic</b>	<b>SDS</b>	<b>0.1%</b>	<b>Yes</b>
	<b>Sodium Deoxycholate</b>	<b>0.5%</b>	<b>Yes</b>

### **Others**

<b>NaCl</b>	<b>1 M</b>	<b>Yes</b>
<b>Glycerol</b>	<b>10%</b>	<b>Yes</b>
<b>EDTA</b>	<b>10 mM</b>	<b>Yes</b>

The “Yes” indicates the reagents can be used in the Lysis buffer for this cartridge up to the indicated concentration. The “No” indicates the reagents cannot be used in the Lysis buffer for this cartridge at the indicated concentration.

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